

QUANTITY	CONTENTS	CODE
30 mL	Cresol Red Indicator	2212-G
2	Test Tubes, 2.5-5-10 mL, plastic, w/caps	0106
1	Cresol Red Octa-Slide 2 Bar, 7.2-8.6	2186-01
1	Octa-Slide 2 Viewer	1101

To order individual reagents or test kit components, use the specified code number.

*Reagent is a potential health hazard. **READ SDS:** [lamotte.com](https://www.lamotte.com)

Emergency information:
Chem-Tel USA 1-800-255-3924
Int'l, call collect, 813-248-0585


Warning! This set contains chemicals that may be harmful if misused. Read cautions on individual containers carefully. Not to be used by children except under adult supervision.




USE OF THE OCTA-SLIDE 2 VIEWER

PROCEDURE


1. Insert Cresol Red Octa-Slide 2 Bar [2186-01] into the Octa-Slide 2 Viewer [1101].

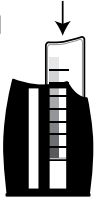


2. Fill a test tube [0106] to the 10 mL line with sample water.




3. Add 10 drops of Cresol Red Indicator [2212].



4. Cap and mix.

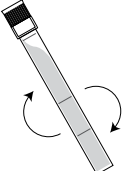

5. Insert test tube into Octa-Slide 2 Viewer [1101]. Hold the Viewer so that non-direct light enters through the back.



6. Match sample color to a color standard. Record as pH.

