

THYMOL BLUE KIT

Code 2112-01 | Octa-Slide 2, pH 8.0-9.4



QUANTITY	CONTENTS	CODE
30 mL	Thymol Blue Indicator	2213-G
2	Test Tubes, 2.5-5-10 mL, plastic, w/caps	0106
1	Thymol Blue Octa-Slide 2 Bar, 8.0-9.4	2187-01
1	Octa-Slide 2 Viewer	1101

To order individual reagents or test kit components, use the specified code number.

*Reagent is a potential health hazard. **READ SDS:** lamotte.com

Emergency information:
Chem-Tel USA 1-800-255-3924
Int'l, call collect, 813-248-0585



Warning! This set contains chemicals that may be harmful if misused. Read cautions on individual containers carefully.
Not to be used by children except under adult supervision.

HOW TO USE THE

OCTA-SLIDE 2 VIEWER

PROCEDURE

1.

Fill a test tube [0106] to the 10 mL line with sample water.

2.

Add 10 drops of Thymol Blue Indicator [2213]

3.

Cap and mix.

4.

Insert Thymol Blue Octa-Slide 2 Bar [2187-01] into the Octa-Slide 2 Viewer [1101]

5.

Insert test tube into Octa-Slide 2 Viewer [1101].

6.

Hold the Viewer so that non-direct light enters through the back. Match sample color to a color standard. Record as pH.